

## 912-0012 **NC-250 IQ/OQ Test Kit**

- P/N: 942-0001** 1 box of **NC-Slide A2** containing 25 glass slides.
- P/N: 071-0015** 1 bottle with 1 mL of fluorescent beads.  
The solution contains multicoloured 10 µm beads,  
Tween-20 (0.02%) and NaN<sub>3</sub> (2 mM ~ 0.01%) in H<sub>2</sub>O.  
CAS no. for Tween-20 is 9005-64-5  
EINECS no. for NaN<sub>3</sub> is 247-852-1

- Application** **NC-250 IQ/OQ Test Kit** is part of the *NucleoCounter*<sup>®</sup> *NC-250*<sup>™</sup> system, and is used for performing installation qualification and operation qualification (IQ/OQ).
- Principle** **NC-250 IQ/OQ Test Kit** contains multicoloured beads and slides that allow testing of the performance of all the channels of the *NucleoCounter*<sup>®</sup> *NC-250*<sup>™</sup>.
- Usage** With respect to the detailed procedure, please refer to the appropriate technical notes and user manuals for the *NucleoCounter*<sup>®</sup> *NC-250*<sup>™</sup> instrument and the *NucleoView NC-250 Software*<sup>™</sup>.
- NC-250 IQ/OQ Test Kit** is for research and development purposes only and is not for diagnostic or therapeutic use.
- Storage** The fluorescent bead solution should be store at 2-7°C. The *NC-Slide A2* should be stored at room temperature < 40°C.
- Stability** The stability of the kit is 12 months if stored under recommended conditions. Expiry dates are shown on the slide box as well as on the fluorescent bead bottle. The **NC-250 IQ/OQ Test Kit** is produced 12 month before the expire date.
- Discard the **NC-250 IQ/OQ Test Kit** after performing the OQ procedure.
- Safety Information** Fluorescent bead solution: In case of contact with eyes, rinse immediately with plenty of water and seek medical advice.  
If a spill is observed perform a clean-up of the area which may have been in contact with the solution.  
  
Use gloves and suitable protective clothing.  
  
Please refer to MSDS regarding further safety information.
- Disposal of Waste** After use the **NC-250 IQ/OQ Test Kit** should be disposed of according to national or local laws and regulations.