

910-3005 Solution 5 - VB-48™•PI•AO**1 ml**

Contents	<p>Solution 5 - VB-48™•PI•AO (1 ml).</p> <p>The solution contains VitaBright -48™ (VB-48™; 400 µg/ml ~ 1.34 mM ~ 0.04%), propidium iodide (PI; 500 µg/ml ~ 0.75 mM ~ 0.05%) and acridine orange (AO; 1.2 µg/ml ~ 3.2 µM ~ 0.00012%) in dimethyl sulfoxide (DMSO).</p> <p>The EINECS no. for PI is 247-081-0.</p> <p>The EINECS no. for AO is 253-353-6.</p> <p>The EINECS no. for DMSO is 200-664-3.</p>
Application	<p>Solution 5 - VB-48™•PI•AO enables detection of changes in cell vitality.</p>
Principle	<p>Solution 5 - VB-48™•PI•AO contains three different dyes: AO stains all nucleated cells, PI stains non-viable cells only, and VB-48™ stains cells in an intensity-dependent manner reflective of their redox state. A high fluorescence intensity of a particular cell indicates that the cell has a high level of reduced thiols. VB-48™ is used to detect mid- to late-stage apoptosis in mammalian cells.</p>
Use	<p>Thaw and mix before use!</p> <p>The solution is for research use only (RUO) and is not for diagnostic or therapeutic uses.</p>
Storage	<p>Store at < 5°C.</p> <p>Protect the solution against moisture and light.</p>
Stability	<p>Unopened microtubes can be stored until the expiry date indicated on the microtube. The solution is produced 15 months prior to the expiry date.</p> <p>We recommend using the solution no later than the expiry date or one (1) month after opening the microtube, whichever comes first.</p>
Safety information	<p>DMSO is an irritating organic compound with long-term effects. DMSO's hazard classification is: Xi.</p> <p>Risk-phrases: R36/37/38 (Irritating to eyes, respiratory system, and skin).</p> <p>Safety-phrases: S26 (In case of contact with eyes, rinse immediately with plenty of water and seek medical advice). S37 (Wear suitable gloves).</p> <p>Please also refer to the safety data sheet (SDS) for safety information.</p>
Disposal	<p>After use, dispose of the bottle following national or local laws and regulations, according to the nature of its contents.</p>